## QUICK START GUIDE

ba77021y01 02/2014

# pH/Cond 3320

a xylem brand

For the most recent version of the manual, please visit  
[www.WTW.com](http://www.WTW.com).

**Copyright**

© Weilheim 2014, WTW GmbH  
Reproduction in whole - or even in part - is prohibited without the  
express written permission of WTW GmbH, Weilheim.  
Printed in Germany.

## pH/Cond 3320 - Contents

EN English .....	4
DE Deutsch .....	6
CS Český .....	8
DA Dansk .....	10
EL Ελληνικά .....	12
ES Español .....	14
FI Suomi .....	16
FR Français .....	18
HU Magyar .....	20
IT Italiano .....	22
JA 日本語 .....	24
KO 한국어 .....	26
NL Nederlands .....	28
NO Norsk .....	30
PL Polski .....	32
PT Português .....	34
RO Română .....	36
RU Русский .....	38
SK Slovenský .....	40
SV Svenska .....	42
TR Türkçe .....	44
ZH 中文 .....	46

## pH/Cond 3320

### Overview

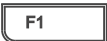
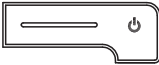
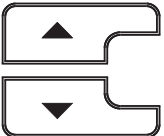
The pH/Cond 3320 meter enables you to perform measurements (pH, conductivity) quickly and reliably.

The pH/Cond 3320 meter provides a maximum degree of operating comfort, reliability and measuring certainty.

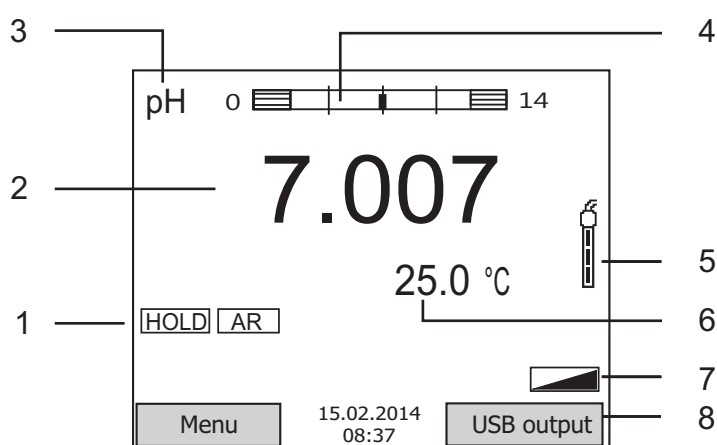
### Keypad

Keys are indicated by brackets <..>.

The key symbol (e.g. <ENTER>) generally indicates a short keystroke (under 2 sec). A long keystroke (approx. 2 sec) is indicated by the underscore behind the key symbol (e.g. <ENTER\_\_>).

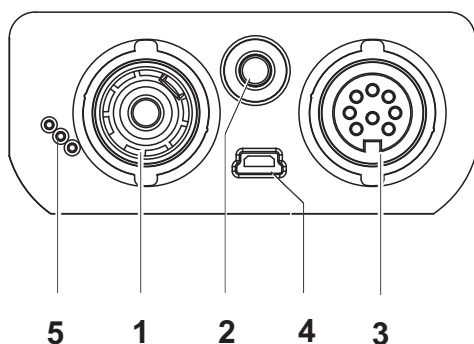
	<F1>: <F1__>:	Situation dependent function (is shown on the display)<F1>/ [Menu]: Opens the menu for measurement settings
	<F2>: <F2__>:	<F1__>/[Menu]: Opens the menu for system settings
	<On/Off>:	Switches the meter on or off
	<M>:	Selects the measured parameter
	<CAL>: <CAL__>:	Calls up the calibration procedure Displays the calibration data
	<STO>: <STO__> :	Saves a measured value manually Opens the menu for the automatic save function
	<RCL>: <RCL__>:	Displays the manually stored measured values Displays the automatically stored measured values
	<▲><▼>:	Menu control, navigation
	<ENTER>: <ENTER__>:	Opens the menu for measurement settings / confirms entries Opens the menu for system settings
	<AR>:	Freezes the measured value (HOLD function) Switches the AutoRead measurement on or off

## Display (example: pH measurement)

1	Status information
2	Measured value (with unit)
3	Measured parameter
4	Continuous measurement control (CMC function)
5	Sensor symbol (calibration evaluation)
6	Measured temperature (with unit)
7	Further status information
8	Softkeys and date/time

## Jack field /connections

1	pH electrode
2	Temperature sensor
3	Conductivity measuring cell
4	USB B interface (device)
5	Service interface

## Initial commissioning

- 1 Insert the supplied batteries.  
Make sure that the poles are in the correct position.
- 2 Switch on the meter with the <On/Off> key.

**Caution**

Follow the safety instructions of the used sensors.  
The operating manuals of the sensors are available on the supplied CD and on the Internet under [www.WTW.com](http://www.WTW.com).

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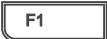
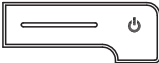
# Xylem |'zīləm|

- 1) The tissue in plants that brings water upward from the roots;
- 2) a leading global water technology company.

We're 12,500 people unified in a common purpose: creating innovative solutions to meet our world's water needs. Developing new technologies that will improve the way water is used, conserved, and re-used in the future is central to our work. We move, treat, analyze, and return water to the environment, and we help people use water efficiently, in their homes, buildings, factories and farms. In more than 150 countries, we have strong, long-standing relationships with customers who know us for our powerful combination of leading product brands and applications expertise, backed by a legacy of innovation.

**For more information on how Xylem can help you, go to [www.xylem.com](http://www.xylem.com)**

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## **Wissenschaftlich-Technische Werkstätten GmbH**