

Geotech Field Equipment Cleaner is a cutting-edge anionic solution formulated to safely decontaminate field and laboratory equipment used for sample collection and analysis. Our FEC solution efficiently eliminates both organic and inorganic contaminants without compromising sample integrity.

## FEATURES

- Complies with EPA Field Equipment Decontamination Guidelines
- Safe, non-corrosive, and non-hazardous formula for worry-free use
- Concentrated formula for efficient cleaning
- Free from 1,4 Dioxane, PFAS, Phosphates
- Leaves no residue after rinsing

## USAGE:

For routine cleaning and decontamination tasks, dilute FEC at a 1:100 ratio with water (1.28 oz per 1 US gallon).

Simply immerse all wetted parts in the diluted solution, dislodge any particulate matter with a brush if necessary, rinse thoroughly with clean tap water, and perform a final rinse with deionized water for optimal cleanliness.

## PERFORMANCE:

Geotech Field Equipment Cleaner is a safe and effective solution that aligns with published environmental field equipment cleaning and decontamination guidelines (LSASDPROC-205-R4). It can be confidently used on all test materials, including metals, glass, and plastics. Please contact Geotech Environmental Equipment, Inc for further information.

## SAFETY WARNING:

- May cause skin irritation.
- Can result in serious eye damage.

## PRECAUTIONARY MEASURES:

- Always wear protective gloves, clothing, eye protection, and face protection when handling.
- If product comes into contact with skin, wash thoroughly with plenty of water.
- In case of eye contact, rinse cautiously with water for several minutes. Remove contact lenses if present and easy to do so, then continue rinsing.
- Immediately contact a poison control center or doctor.
- Remove contaminated clothing and wash before reuse.

Prioritizing safety is paramount when handling this product. Please adhere to the precautionary statements for your well-being and protection.



CALL GEOTECH TODAY

**(800) 833-7958**

Geotech Environmental Equipment, Inc.

2650 East 40th Avenue Denver, Colorado 80205

(303) 320-4764 FAX (303) 322 7242

email: sales@geotechenv.com website: www.geotechenv.com